



## Microlucida®

### Release Notes

## Version 2026.1.1 (February 2026)

### NEW FEATURES AND ENHANCEMENTS

- New Background Subtraction and Flat Field Correction image filters added to Slide Scanning and image compilation
- New Blending function in the Slide Scanning workflow
- Automatic stitching of overlapping fields of view in the Montage tool is now available
- AutoMove will now reposition focus when Continuous tracing is being used
- Image file paths are displayed in the Image Organizer

### ISSUE RESOLVED

- Fixed crash in Slide Scan workflow sample scan
- Optimized settings for BSI Prime camera

## Version 2025.1.3 (July 2025)

### ISSUE RESOLVED

- Fixed licensing issue that would prevent the program from running
- Fixed a crash in slide scanning choosing to acquire stack at a time while alternating channel order

## Version 2025.1.2 (May 2025)

### NEW FEATURES AND ENHANCEMENTS

- Support added for the new Lumina 9 camera
- Improvements made to image acquisition with the Hamamatsu camera

### ISSUES RESOLVED

- Fixed an issue that could occur when using the Apotome during a slide scan image acquisition
- The storage estimate needed for stitching during a slide scan was fixed

## Version 2024.1.3 (November 2024)

### LICENSING UPDATE

- Update to offline software licensing. MicroLucida software requires reactivation on systems that do not have access to the internet. [Instructions for offline software activation](#) are included in the MicroLucida user guide.

### ISSUES RESOLVED

- Fixed a problem with Live Image not working when the program is started
- Fixed an issue with the Live Zoom button on the Acquire ribbon