



NeuroInfo® software

Release Notes

Version 2025.1.1 (November 2025)

NEW FEATURES AND ENHANCEMENTS

- New Map to Intermediate option is available after completing Section Registration using Atlas-constrained registration to create a brain image volume registration that can be further refined for better precision using the Register Volume tools
- New Nonlinear Registration options added provide better results and more control of the parameters used
- Calibration and transform information from image registration are now stored with the data file
- Additional tools and information in the Calibration panel:
 - View information about installed atlases
 - Atlas names now include the species, version, and section spacing
 - Activate an updated or different brain atlas to modify the transforms from a previous registration
 - New warning if the atlas and/or spacing is not compatible with the experimental images
- Improvements to the Tone Mapping filter
- New marker density visualization surface option added to the 3D Visualization window
- Large image and data file handling are now faster and more robust
- Improvements and updates to the Movie mode interface have been implemented
- The Gerfen Nissl Mouse Atlas has been updated to version 2

ISSUES RESOLVED

- Resolved resolution issue during the Serial Section Assembler setup

Version 2024.1.3 (December 2024)

LICENSING UPDATE

- Update to offline software licensing. NeuroInfo software requires reactivation on systems that do not have access to the internet. [Instructions for offline software activation](#) are included in the NeuroInfo user guide.

NEW FEATURES AND ENHANCEMENTS

- Automove will now reposition focus when Continuous tracing is being used
- Improvements to synchronization of contours delineated in 2D and 3D
- Image file paths are displayed in the Image Organizer
- During a Movie, you can now change image-display settings for image slice and partial projections
- New ability to display the 3D scale bar while recording a Movie
- Other improvements and updates to the Movie mode interface have been implemented

ISSUES RESOLVED

- Resolved an issue exiting quad view
- An image will close properly if one channel was selected prior to closing all images