



Neurolucida® software

Release Notes

Version 2021.1.1 (5/25/2021)

- Add ability to rotate individual points on a contour/tree
- Added the ability to end tree tracing early, with a right click option. Added the ability to resume tracing by clicking on a node
- Support for new Lumenera Lt-3200 CMOS color camera
- Support for Zeiss MTB version 2.17.0.15
- Support for Nikon RDK 1.4.0.115

ISSUES RESOLVED

- If the Hamamatsu camera becomes unresponsive during an acquire the software will now gracefully end the acquisition enabling the post acquire command sequences to run including closing the shutter, if configured
- Fixed a bug in slide scan acquires that would cause a crash if you entered a non-standard file extension (not jpeg 2000) when entering the image file name in the workflow
- When doing slide scans, write out settings to a text file (MyImg_AcquireSettings.txt) so we can see what they were at the time of the acquire
- Fixed a bug in the slide scanning workflow where users could get stuck with a bad step size value in 2D scans