



# Stereo Investigator®

Release Notes

## Version 2022.3.1 (December 2022)

### NEW FEATURES AND ENHANCEMENTS

- New *Batch Pipeline* enables easy-to-use batch processing for image filtering of neuronal structures
- New *Pipelines* ribbon now houses features previously located in the *Automatic* section of the *Trace* Ribbon
- Improved & expanded image support
- New automatic, smart image-optimization preference for low-signal images available
- The subvolume tool is now restricted by the bounds of the image
- Image montage can now assemble images of varying x/y sizes

### ISSUES RESOLVED

Numerous bug fixes throughout the software. These include:

- Fixed options for saving the current plane
- Fixed a problem with software becoming unresponsive when acquiring images with Microfire cameras
- Fixed crashing associated with acquiring images with Apotome
- Fixed crashing when the modal version of the exposure dialog bar was selected during a background correction step in the Slide Scanning workflow
- Fixed a problem with the save location being ignored when post-compiling a Slide Scan with preview and stitching options selected.
- Fixed a crash that occurred rarely during slide scanning workflows
- Fixed a crash that occurred in specific conditions when the stage is moved with the Image Organizer open.