



# Stereo Investigator Whole Slide Edition

Release Notes

## Version 2020.1.1 (2/20/2020)

### NEW FEATURES AND ENHANCEMENTS

- Added controls in the export tracing dialog for exporting images at a specific resolution
- Added an edit option to 'Select everything of this contour type'
- Added a new, fast cursor option in preferences which addresses potential cursor delay issues
- Shift click will now erase markers in Point Sampled Intercept probe
- New montage option to 'Order and align using image name'

### ISSUES RESOLVED

- Fixed image export issue that exported the background image at a low resolution.
- Fixed issue with Physical Fractionator when re-opening a previously completed run:
  - First problem caused a crash
  - Second problem prevented the user from moving past the step
- Fixed issue where rotating jpeg2000 images could result in images that were too large or in the wrong location.
  - If running Physical Fractionator, this can result in the counting frame not landing on the correct part of the rotated image
- Fixed issue with Physical Fractionator on a 4K monitor
- Fixed crash that occurred when cancelling the tool panel "Close All" button with Image Adjustment window
- Fixed issue where the lines in the Merz probe were too thin on a 4k monitor
- Fixed issue with transparency slider in preferences
- Fixed issue when loading RGB nd2 images
- Fixed issue where scaling is over-ridden when images are opened in grid layout
- Fixed issue that 'Show current section' was not respected in the 3D visualization window
- Fixed crash that occurred when changing the tracing options with the circle contour